

dH2O rinse (10 min)			
dH2O rinse (10 min)			
dH2O rinse (10 min)			
Uranyl acetate stain 1hr			
dH2O rinse (10 min)			
dH2O rinse (10 min)			
dH2O rinse (10 min)			
50% EtOH (10-20 min)			
50% EtOH (10-20 min)			
70% EtOH (10-20 min)			
70% EtOH (10-20 min)			
80% EtOH (10-20 min)			
80% EtOH (10-20 min)			
90% EtOH (10-20 min)			
90% EtOH (10-20 min)			
100% EtOH (10-20 min)			
100% EtOH (10-20 min)			
100% EtOH (10-20 min)			
Resin/Propylene Oxide (1:1) (1h-O/N)			
Embedding Resin (1-2 h)			
Embedding Resin (1-2 h)			
Resin curing (60°C, O/N)			